
Visit our website I Visit our blog I Like us on facebook

■ instrument choice.com.au[™]

July 2013

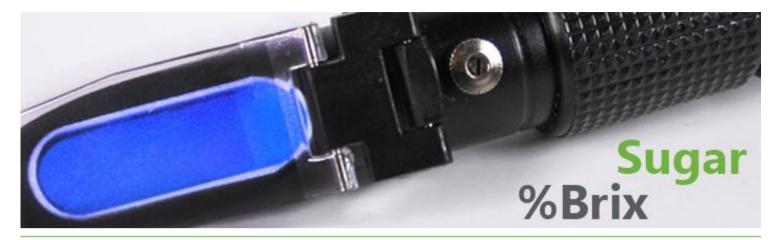

July 2013 Newsletter - Experimenting with Refractometers and Honey.

We have some pretty exciting news to share this month. Instrument Choice (Synotronics Pty Ltd) has been announced as a finalist in this year's <u>Telstra Australian Business Awards</u>. We were lucky enough to be finalist last year and this year we are hoping for the win. We are up against some very good business' so it won't be easy. You can check out our competition on the <u>Australian Business Awards website</u>.

We will keep you up dated of how we go, and we will be facebooking live from the event, you can follow us <u>here</u>. Win or lose we are all really looking forward to the Gala Dinner, it is a great night and a chance for everyone to let their hair down after another big year.

Over the past couple of months we have noticed an increase in the number of <u>refractometer</u> enquiries we have been receiving. This month Daniel puts a <u>Honey refractometer</u> through its paces, testing a range of supermarket bought honeys. We have been madly adding new products to our range over the past couple of weeks so Tim will be introducing you to some of our new standout lines.

Remember; if you have any questions regarding tests you need to conduct or specific instrumentation please give us a call.

Until next month... Tyson Grubb Product of the month:

HI96801 - Digital

Refractometer for Sugar



A Digital refractometer ideal for the analysis of fruits, energy drinks, puddings, soy milk juices, jam, marmalade, honey, soups, jelly, tofu and condiments.

The dual-level LCD displays measurement and temperature readings simultaneously along with "Automatic Temperature Compensation" for accurate measurements.



Sweets for my sweet and sugar for my honey.

This month we will be testing different types of honey with a refractometer to determine their water content. Honey with a water content greater than 20% is at a higher risk of fermentation. Fermentation causes a reduction in honey quality and adversely affects the taste. Read on to find out what we did!

Free Express Post in Australia. Items marked bulky have free standard post.






Equipment used:

- IC-MR90ATC Honey refractometer
- 6 types of honey
- Beaker
- Demineralised water
- Teaspoons
- Olive oil

Lint free cloth Ceramic dinner plate

Method:

Each of the honey containers were opened and a small sample was placed on a ceramic plate ready for measurement. The IC-MR90ATC refractometer was calibrated to 72% Brix with olive oil (72% is the known Brix level of olive oil) prior to the initial test to ensure accuracy. After calibration, each of the samples were measured for their water content and this was recorded. We also recorded the Brix and Baume readings for reference. The refractometer was cleaned with a soft cloth and demineralised water between each measurement. We also recorded the temperature and humidity of the room prior to each reading for each of the samples for reference using a hygrometer.

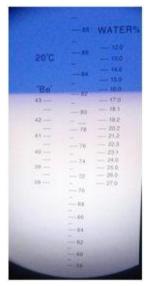

View through refractometer eye piece for Blue gum honey

Samples tested:

We tested 6 types of honey to get a good gauge on the types of moisture content that are acceptable in packaged honey. The samples included Blue gum, Kangaroo Island, Leatherwood, Creamed, Home brand, and Fast food.

Results:

Sample	Blue Gum	Kangaroo Island	Leatherwood	Creamed	Home brand	Fast food sachet
Moisture Content %	16.2	17.5	16.8	N/A	17	16.5
Brix %	82.1	80.9	81.8	N/A	81	81.9
Baume	Off scale	42.7	Off scale	N/A	42.8	Off scale
Temperature (°C)	22.6	22.6	22.6	22.6	22.6	22.6

Humidity (%) 41.1 42 40.7 40.7 41.2 40.7

Discussions and conclusions:

All of the samples we tested were well below 20% moisture content and as such are at less risk of fermentation. The honey with the lowest moisture content was the Blue Gum honey at 16.2%, whilst the highest was the Kangaroo Island Honey at 17.5%. It is interesting to note that all of the honeys are within 1.3% of each other. We were unable to obtain a result from the creamed honey as it kept giving erroneous values. This was due to the fact it is an opaque liquid that wouldn't allow the light to pass through like the other honeys and therefore would not give a correct moisture or Brix measurement.

Honey with a high moisture content is at risk of fermenting due to the presence of yeast species within the honey. At low moisture contents these yeasts cannot grow as the sugar depletes the water from the yeast cells. However, as the moisture content increases the yeast cells no longer remain dormant and can begin to grow. As the yeast cells multiply they produce alcohol. The produced alcohols can then break down and turn into acetic acid and water. This gives honey a sour and off taste. It may also result in a change in texture of the honey, small bubbles within the honey or foaming on the surface.

It also interesting to note that honey is a hygroscopic liquid, meaning it attracts moisture from the air. Therefore in a very dry, warm atmosphere the moisture content will actually decrease if it is exposed to this environment and vice versa if it is exposed to a cold, moist environment. Therefore keeping your honey stored in an air tight container will lessen the risk of fermentation as it cannot be affected by moisture in the air.

Hopefully you have found this month's newsletter enjoyable, if you have any questions please <u>contact</u> one of our scientists and they will be only too happy to assist you.



What's new this month?

Every month we are adding new exciting products to the website, here are the latest for July.

Casella dBadge personal sound exposure meters

Horiba range of water quality meters

Dickson digital chart recorders

This month has seen the addition of the <u>Casella</u> dBadge personal sound exposure meters to our website. The dBadge dosimeters are compact and rugged and are light enough to be worn on the shoulder. The standard <u>dBadge measurement kit</u> contains everything you need to start measuring personal noise exposure in the workplace and will allow you to view cumulative noise dose data and view time history profiles using the software provided.

The dBadge Plus has the same features as the standard dBadge but also allows for the simultaneous measurement of LCeq and LAeq which can be used for selection of suitable hearing protection. Intrinsically safe kits of both the dBadge and dBadge Plus are also available and it is worth keeping in mind that every Casella sound level meter, logger, dosimeter and calibrator comes with a signed and dated calibration certificate from the manufacturer.



We have also added the Horiba range of water quality meters to our website. Horiba manufacture a wide range of high quality water quality meters from compact single parameter meters through to multi-parameter laboratory and field instruments.

The compact LAQUA range of single parameter testers allow you to accurately take measurements of pH, conductivity and salt as well as sodium, potassium, nitrate and calcium ions from a single drop of solution. The B-711 has already been popular with customers measuring the pH of plant sap. For customers looking to measure dissolved oxygen or Biochemical Oxygen Demand (BOD), the OM-51-L1 is a great option as the electrode has a standard taper designed for use with

standard BOD bottles.

For those of you that are interested measuring temperature and humidity and considering in moving from a paper chart recorder over to a digital version, the Dickson digital chart recorders could be just the ticket. We currently have the FH525 and the FH635 models on the website. The FH525 has a temperature and humidity sensor on a 3m cable and the FH635 has a temperature and humidity sensor which is built into the unit.

They both have a 104 x 140mm touchscreen display as well as audio and visual alarms. The FH635 also features 2 SPST 24VDC 0.5A NO/NC relays. The touch screen display has full graphing capabilities so you can quickly view temperature and humidity history. Optional software is also available if you want to download the data to your PC.



Thank you

from everyone at Instrument Choice - stay tuned for next month's issue.



Contact us.

Our experts are happy to help and discuss your project. Call 1300 737 871 or write an email to customer-service@instrumentchoice.com.au















