## HALO 2

# Wireless Refillable pH Tester for Cosmetic Creams

with built-in specialized electrode



# INSTRUCTION MANUAL



## Dear Customer,

Thank you for choosing a Hanna Instruments product.

Please read this instruction manual carefully before using this instrument.

This manual will provide you with the necessary information for correct use of this instrument, as well as a precise idea of its versatility.

If you need additional technical information, do not hesitate to e-mail us at tech@hannainst.com or view our contact list at www.hannainst.com.

## **TABLE OF CONTENTS**

1. Preliminary Examination	4
2. Specifications	4
3. General Description & Intended Use	5
Operating Modes	
Probe Features	5
4. Functional Description & LCD Display	5
5. General Operations	6
Turning the Tester On & Off	6
Battery Replacement	6
6. Setup	
Temperature Unit	
Auto-Off Interval	
Calibration Points	
Buffer Set	
pH Resolution	
Bluetooth Pairing	
Date & Time	
7. Bluetooth	
8. Hanna Lab Application	
Tag a Measurement	
9. Calibration	
Preparation & Guidelines	7
Procedure	
Stand-Alone Tester (Up to Three-Point Calibration)	
With Hanna Lab App (Up to Four-Point Calibration)	
10. Care & Maintenance	
Refilling the Electrode	
Storage	
11. Warning & Error Messages	
12. Abbreviations	
13. Accessories	
Recommendations for Users	
Warranty	
Certification	11

## 1. PRELIMINARY EXAMINATION

Remove the tester and accessories from packaging and examine them carefully. For further assistance, please contact your local Hanna Instruments office or email us at tech@hannainst.com.

Each H19810432 is supplied with a starter kit consisting of:

- pH 4.01 buffer solution, 20 mL sachet (2 pcs.)
- pH 7.01 buffer solution, 20 mL sachet (2 pcs.)
- Electrode cleaning solution, 20 mL sachet (2 pcs.)
- Electrode storage solution, 13 mL dropper bottle (1 pc.)
- Electrolyte refill solution, 30 mL
- Pipette
- 3V Lithium battery CR2032
- Instrument quality certificate and Instruction manual

**Note:** Save all packing material until you are sure that the tester works correctly. Any damaged or defective item must be returned in its original packing material with the supplied accessories.

## 2. SPECIFICATIONS

рН	0.00 to 12.00 pH	
mV *	pH/mV conversion	
Temperature**	−5.0 to 70.0 °C (23.0 to 158.0 °F)	
рН	0.01 or 0.1 pH	
mV *	0.1 or 1 mV	
Temperature	0.1 °C; 0.1 °F	
pН	$\pm 0.02  \mathrm{pH}$	
Temperature	$\pm 0.5$ °C; $\pm 0.9$ °F	
Automatic buffer	ts or four points * recognition with Standard buffers t *, 4.01, 7.01, 10.01) or NIST (pH 1.68 *, 4.01, 6.86, 9.18)	
Automatic (ATC)	or Manual (MTC) *	
Body material Glass Junction Reference cell Electrolyte Tip / Shape Outer diameter Length	Glass Low Temperature (LT) Triple ceramic Double, Ag/AgCl 3.5M KCl (refillable) Conic 12 mm (0.47") 120 mm (4.7")	
3V Lithium — CR	2032	
Approximately 1000 hours (500 hours with Bluetooth enabled)		
0 to 50 °C (32 to 122 °F)		
IP65 ingress pro	tection	
51 v 206 v 21 m	nm (2.0 x 8.1 x 0.8") / 60 g (2.1 oz.)	
	mV * Temperature** pH mV * Temperature pH Temperature Up to three poin Automatic buffer Hanna (pH 1.68 Automatic (ATC) Body material Glass Junction Reference cell Electrolyte Tip / Shape Outer diameter Length 3V Lithium — CR Approximately 1 0 to 50 °C (32 t IP65 ingress pro	

<sup>\*</sup> Available with Hanna Lab App

<sup>\*\*</sup> Measuring outside the recommended operating temperature range may damage the gel electrolyte and void product warranty.

Note: The tester can display measurements from -2.00 to 16.00 pH. Measurements outside of the pH range will flash. In this case assess the integrity of the tester and the type of measured sample.

## 3. GENERAL DESCRIPTION & INTENDED USE

HI9810432 is a professional wireless pH tester, part of Hanna Instruments HALO2 family.

- The integrated Bluetooth module allows the tester to be connected to a compatible smart device with the Hanna Lab App.
- The tester has a compact, waterproof casing, and automatic pH calibration at up to three points, or four points when
  used with the Hanna Lab App.
- Automatically compensated temperature readings are displayed on a large LCD.
- Accurate and easy to use, the tester is ideal for measurements in samples that would be a challenge for standard design pH electrodes.

## **Operating Modes**

The H19810432 can be used as a stand-alone pH tester or connected to the Hanna Lab App.

The Hanna Lab App turns a compatible smart device into a full-featured pH meter. Features include: electrode condition, GLP with time-stamp, live-readings, mV resolution, manual temperature compensation, stability criteria, calibration reminder, pH (mV) and temperature alarms, tester ID, and data sharing.

## **Probe Features**

The HI9810432 features a non-porous, glass body that is easy to clean and withstands harsh chemicals.


The conical glass tip allows for penetration into semisolids and emulsions.


The **triple ceramic junction** allows a higher flow rate of electrolyte from the reference cell into the solution. This high flow rate provides faster electrode response and a more stable measurement in viscous solutions or samples of low conductivity.

Built-in **temperature sensor** at the tip of the pH electrode allows for rapid determination of the sample temperature and a high-accuracy temperature reading.


## 4. FUNCTIONAL DESCRIPTION & LCD DISPLAY

## Front & Back view

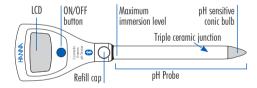

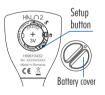

## LCD display

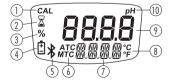

- 1 Calibration indicator
- 2 Stability indicator
- 3 Battery percentage indicator
- 4 Battery icon
- 5 Bluetooth icon
- 6 Temperature compensation
- 7 Second LCD line
  - 8 Temperature unit
- 9 First LCD line
- 10 Measurement unit

## 5. GENERAL OPERATIONS

## Turning the Tester On & Off

- Turn the tester face down on a clean, dry surface. Rotate the battery cover on the back of the tester counterclockwise and set it aside. Remove the battery insulation film.
- Press the ON/OFF button to turn the tester on or off. Initialization screen displays all LCD segments, followed by the
  battery percentage. The tester enters measurement mode. Before the tester turns off, "BFF FWR" is displayed briefly.
- Press and hold the ON/OFF button to turn the tester off when connected to Bluetooth.

## **Battery Replacement**

- 1. Turn off the tester. Turn the tester face down and rotate the battery cover counterclockwise.
- 2. Set the battery cover aside. Press the metallic pin to push old battery out.
- 3. Place the new battery with positive (+) sign facing out.
- 4. Reset the date and time in setup, or connect to the Hanna Lab App to update it automatically.
- Align the mark on the cover with the open lock icon (ar) on the case. Rotate the cover clockwise until the mark on the cover aligns with the closed lock icon (a).

Note: Only use specified battery type. Dispose of old battery in accordance with local regulations.



## 6. SETUP

The Setup button is located inside the battery compartment. After Setup configuration, replace the cover.

## Setup Menu Navigation

- Press the Setup button to enter Setup mode and navigate menu items.
- To exit Setup mode, press the Setup button after "SEL TIME" option is displayed.
- Press the ON/OFF button to configure menu item options.

## **Temperature Unit**

Option: °C or °F
Press ON/OFF button to select desired

temperature unit.

## Auto-Off Interval

Option: 8, 60 min., or "----" (disabled)
Press ON/OFF button to select desired interval.
To save battery life, after the selected auto-off

interval has elapsed, the tester will automatically turn off.



op

## **Calibration Points**

Option: 2P or 3P

Press ON/OFF button to select between up to two- or up to three-point calibration.

## **Buffer Set**

Option: 7.01 pH (Hanna) or 6.86 pH (NIST)
Press ON/OFF button to select the calibration
buffer set (Hanna or NIST).

## pH Resolution

**Option:** 0.01 pH or 0.1 pH

Press the ON/OFF button to select resolution.

## **Bluetooth Wireless Mode**

Option: On, PAIr, or OFF

Press ON/OFF button to select Bluetooth

option at start-up.



## **Bluetooth Pairing**

Option: dEL PAIr

Press ON/OFF button to delete saved paired device.

## Date & Time

Option: SEt TIME

Press ON/OFF button to set the date and time.
Option: YEAR. MO. DAY. HOUR. and MIN

NE utton to set the date and time









Use the Setup button to select the option and press the ON/OFF button to change the selected option.

## 7. BLUETOOTH

With "PRI r BLL!" or "On BLL!" selected in Setup, the Bluetooth icon (\*) will blink for up to 45 seconds, indicating the tester is in discoverable mode. Once connected, the icon will stop blinking. If not connected, the icon is not displayed.

- Select " ${\it D}_{\it n}$   ${\it BLU}$ ", to enable Bluetooth without bonding.
- Select "PRI = 31,11", to enable Bluetooth with bonding. A 6-digit bonding pin is displayed the first time the tester and smart device are paired. Once the devices are paired, the bonding pin is not required when reconnected.
- Select "OFF BLU", to disable Bluetooth.
- Select "AFL PAIR", to delete all paired devices. If PAIr BLU is enabled, a bonding pin will need to be re-entered.

## 8. HANNA LAB APPLICATION

- The Hanna Lab App is available on the App Store® and on Google Play.
- Consult the Help section of the application for information on calibration, measurement, data logging, and sharing.
- When the tester is in discoverable mode, it will appear in the list of "Available Devices".
- Within the application, tap "Connect" to pair the tester with the device. All readings are transmitted directly to the application.

## Tag a Measurement

Once connected to the application, the ON/OFF button can be pressed to tag the current reading.

- Press the ON/OFF button from measurement mode. The display shows "SEL TRG", followed by "- TRG".
- The reading on the application will flash green and the note icon ((C)) will be displayed. Tap the note icon ((C)) to add an apportation





## 9. CALIBRATION

## **Preparation & Guidelines**

- Unscrew the top collar to vent the cap before removing the probe from the plastic storage tube. Save the cap for probe storage.
- 2. Rinse off any storage solution or salts that may be on the body.
- Verify there is solution inside the pH bulb by shaking the probe down to restore continuity as the solution may have moved up the stem during shipping.
- 4. For best results, use a rinse beaker and a separate calibration beaker for each buffer. Discard rinsing buffers after use.

## Procedure

For most applications it is recommended to start with pH 7.01 buffer (or pH 6.86).

To restore factory defaults, press and hold the ON/OFF button. "**ERL**  $\square \square R$ " is displayed.

Note: It is recommended to calibrate the electrode with buffers at the temperature it will be used at.

App Store is a trademark of Apple, Inc.

## Stand-Alone Tester (Up to Three-Point Calibration)

- 1. Rinse the electrode tip with purified water and blot dry. Then rinse with the buffer being used for calibration.
- 2. Press and hold ON/OFF button until "ERL METTE" is displayed.
- 3. When "7.0 ! USE" or "6.86 USE" is displayed with "CAL" tag blinking, place the tip of the electrode in the correct buffer.



- 4. When the buffer has been recognized, "REE" is displayed.
  - "WATTT" is displayed with the stability indicator (\$\overline{\mathbb{Z}}\) blinking until the reading is stable. Wait until the measurement is stored and the stability indicator disappears.



To save a one-point calibration and return to measurement mode, press the ON/OFF button.
 "ERL SAVE" is displayed briefly.

## With 2P option selected

- Place the tip of the electrode in the second buffer rinse beaker, then in the second calibration buffer. Wait until the measurement is stored and the stability indicator disappears.
- After the second point has been stored, "EAL SAVE" is displayed briefly and the tester will return to measurement mode automatically.

## With 3P option selected

- Place the tip of the electrode in the second buffer rinse beaker, then in the second calibration buffer. Wait until the measurement is stored and the stability indicator disappears.
- Place the tip of the electrode in the third buffer rinse beaker, then in the third calibration buffer. Wait until the measurement is stored and the stability indicator disappears.
- After the third point has been stored, "ERL SAVE" is displayed briefly and the tester will return to measurement mode automatically.

## With Hanna Lab App (Up to Four-Point Calibration)

Connect the tester to the Hanna Lab App and follow the calibration procedure.

See the App Help section for the calibration procedure.

## 10. CARE & MAINTENANCE

- Never immerse the tester over the maximum immersion level.
- Fresh buffers should be used for each calibration. Once the sachets are opened the buffer value can change over time.
- If the electrode is slow or sluggish, soak it in cleaning solution for 20 minutes. Rinse with water and hydrate the electrode
  in storage solution for a minimum of 30 minutes before calibrating.
- If measurements are taken successively, rinse the probe thoroughly in distilled or deionized water to eliminate cross-contamination between measurements.

## Refilling the Electrode

- If the filling solution (electrolyte) is more than 1 cm (½") below the fill hole, add H17082 3.5M KCl electrolyte fill
  solution for double junction electrodes.
- Unscrew the fill hole cover during measurements so the liquid reference junction maintains an outward flow of electrolyte.

## Storage

To ensure a quick response, the glass tip and junction should be kept hydrated.

 When not in use, add a few drops of storage solution to the storage cap. If storage solution is not available, pH 4.01 or pH 7.01 buffer can be used.

Note: Do not store the electrode in distilled or deionized water.

2. Place the probe tip into the cap and gently tighten the collar.

## 11. WARNING & ERROR MESSAGES



"---- WRNG" displayed during calibration.
Invalid buffer.
Check the buffer value and use fresh buffer



Measured value displayed blinking.

Measured value is out of electrode range.

Clean the electrode to improve condition.



Battery icon (由) displayed blinking. Battery is below 10 %, replace the battery.



"BAtt DEAD" and tester turns off. Depleted battery, replace the battery.

## 12. ABBREVIATIONS

ATC Automatic Temperature Compensation

GLP Good Laboratory Practice

MTC Manual Temperature Compensation

NIST National Institute of Standards and Technology

## 13. ACCESSORIES

Ordering Information	Product Description
pH Buffer Solution	1 rodest postriprior
HI50016-02	pH 1.68 buffer solution, 20 mL sachet (25 pcs.)
HI70004P	pH 4.01 buffer solution, 20 mL sachet (25 pcs.)
HI70006P	pH 6.86 buffer solution, 20 mL sachet (25 pcs.)
HI70007P	pH 7.01 buffer solution, 20 mL sachet (25 pcs.)
HI70009P	pH 9.18 buffer solution, 20 mL sachet (25 pcs.)
HI70010P	pH 10.01 buffer solution, 20 mL sachet (25 pcs.)
HI77400P	pH 4.01 & 7.01 buffer solution, 20 mL sachet (10 pcs., 5 each)
HI770710P	pH 10.01 & 7.01 buffer solution, 20 mL sachet (10 pcs., 5 each)
Electrode Cleaning Solu	
HI700601P	Electrode cleaning solution for general use, 20 mL sachet (25 pcs.)
HI7073L	Protein cleaning solution, 500 mL
HI7074L	Inorganic cleaning solution, 500 mL
HI7077L	Oil & Fat cleaning solution, 500 mL
Electrode Storage Solut	•
HI70300L	Electrode storage solution, 500 mL
HI70300M	Electrode storage solution, 230 mL
HI70300S	Electrode storage solution, 30 mL dropper
HI9072	Electrode storage solution, 13 mL dropper
Electrode Fill Solution 8	Cother Accessories
HI7082	3.5M KCl electrolyte fill solution for double junction electrodes (4 x 30 mL)
HI740155P	Electrode refilling pipette (20 pcs.)

## RECOMMENDATIONS FOR USERS

Before using this product, make sure it is entirely suitable for your specific application and for the environment in which it is used. Any variation introduced by the user to the supplied equipment may degrade the tester's performance. For your and the tester's safety do not use or store it in hazardous environments.

## WARRANTY

H19810432 is warranted for a period of one year against defects in workmanship and materials when used for its intended purpose and maintained according to instructions. This warranty is limited to repair or replacement free of charge. Damage due to accidents, misuse, tampering, or lack of prescribed maintenance is not covered. If service is required, contact your local Hanna Instruments office. If under warranty, report the model number, date of purchase, serial number, and the nature of the problem. If the repair is not covered by the warranty, you will be notified of the charges incurred. If the tester is to be returned to Hanna Instruments, first obtain a Returned Goods Authorization (RGA) number from the Technical Service department and then send it with shipping costs prepaid. When shipping any product, make sure it is shipped back in compliance with shipping regulations, thoroughly cleaned and free of any chemicals, and is properly packaged for complete protection.

10 MAN9810432 08/21

## **CERTIFICATION**

All Hanna Instruments conform to the CE European Directives.

Disposal of Electrical and Electronic Equipment. The product should not be treated as household waste. Instead, hand it over to the appropriate collection point for the recycling of electrical and electronic equipment. which will conserve natural resources.



**RoHS** compliant



**Disposal of waste batteries.** This product contains batteries, do not dispose of them with other household waste. Hand them over to the appropriate collection point for recycling. Ensuring proper product and battery disposal prevents potential negative consequences for the environment and human health. For more information, contact your city, your local household waste disposal service, or the place of purchase.

Regulatory Notices for the stand-alone, Bluetooth, low-energy module.

United States (FCC) FCC ID: 2AA9B04. This device complies with FCC Rules. Part 15 Subpart C "Intentional Radiators" and Subpart B. Chapter 815, 105. This equipment has been tested and found to comply with the limits for a Class A digital device, pursuant to part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference when the equipment is operated in a commercial environment. This equipment generates, uses, and can radiate radio frequency energy and, if not installed and used in accordance with the instruction manual, may cause harmful interference to radio communications. Operation of this equipment in a residential area is likely to cause harmful interference in which case, users are required to correct the interference at their own expense.

Canada (ISED) IC: 12208A-04. This device complies with Industry Canada license exempt RSS standard(s). Operation is subject to the following two conditions: (1) this device may not cause interference, and (2) this device must accept any interference, including interference that may cause undesired operation of the device. Le présent appareil est conforme aux CNR d'Industrie Canada applicables aux appareils radio exempts de licence. L'exploitation est autorisée aux deux conditions suivantes: (1) l'appareil ne doit pas produire de brouillage, et (2) l'utilisateur de l'appareil doit accepter tout brouillage radioélectrique subi, même si le brouillage est susceptible d'en compromettre le fonctionnement.

Australia / New Zealand (RCM) BMD-300 complies with the AS/NZS 4268:2017.

Japan (MIC) (R210-106799

South Korea (KCC) R-CRM-Rgd-BMD-300

Brazil (ANATEL): Contains ANATEL approved module # 00820-21-05903.

Mexico (IFETEL): Este equipo contiene el módulo con IFT #: NYCE/CT/0146/17/TS.

## World Headquarters

Hanna Instruments Inc. Highland Industrial Park 584 Park East Drive Woonsocket, RI 02895 USA www.hannainst.com



MAN9810432

Printed in ROMANIA