**i-DT61XWE Indicator
Ethernet Option
Instruction Manual**

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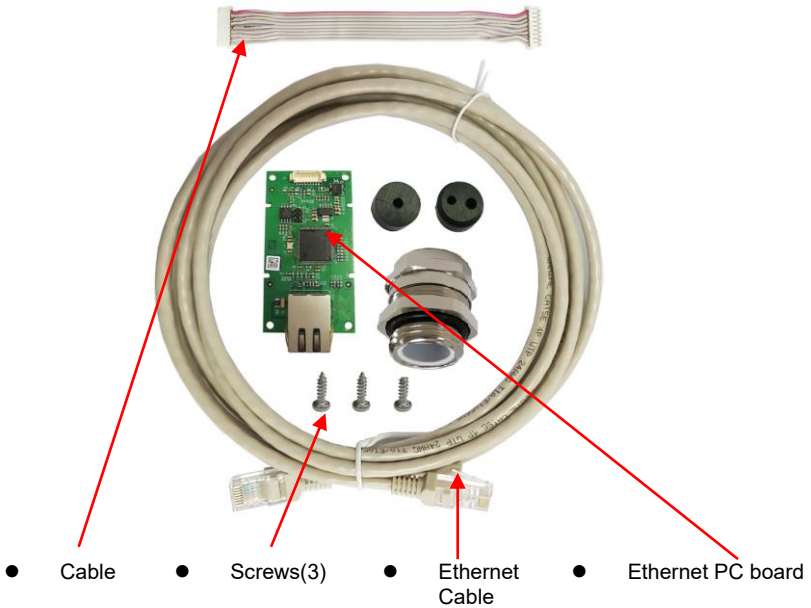
INTRODUCTION

This Ethernet interface kit is for use with the Defender® 6000 i-DT61XWE indicators.

CAUTION: THE INSTALLATION AND CONNECTIONS SHOULD ONLY BE DONE BY AUTHORIZED AND QUALIFIED PERSONNEL, SUCH AS AN ELECTRICIAN.


KIT CONTENTS


INTERFACE INSTALLATION

Install the Ethernet Interface kit to the indicator as the following.

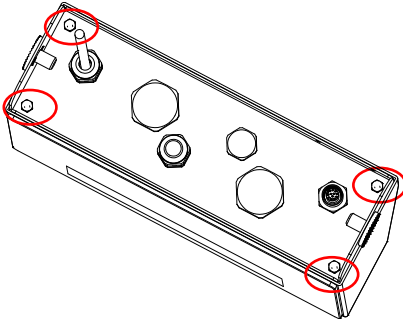
1. Remove the four hex head bolts.

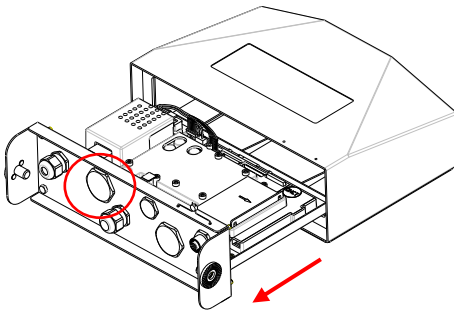
2. Pull out the bottom of the indicator, and remove the following circled M25 cover

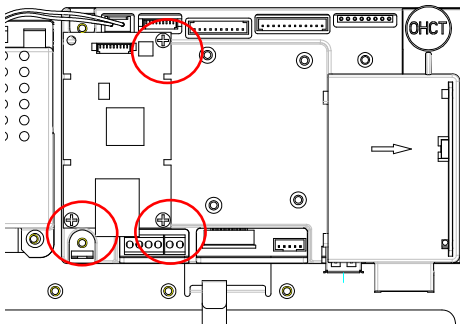
3. Install the Ethernet board. Secure it with 3 screws.

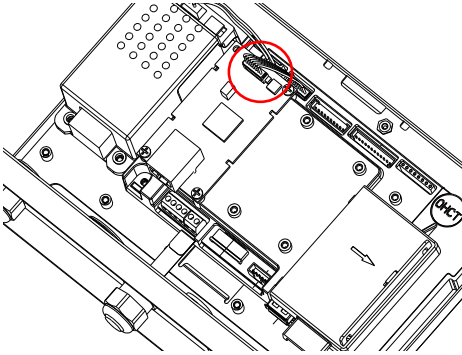
4. Connect the cable to the main board connector.

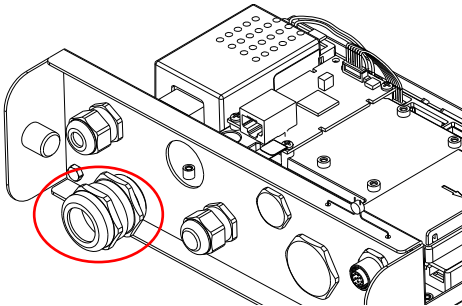
5. Insert the M25 cable gland into the rear housing.

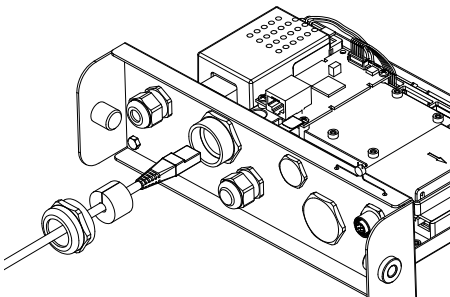6. Disassemble the cable gland by loosening the cap nut and removing the rubber seal. Pass the Ethernet cable through the cap nut, then spread the rubber seal and place it over the cable. After that, pass the cable through the cable gland.

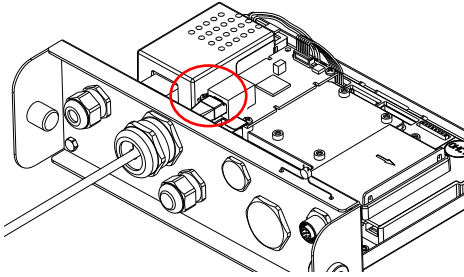7. Connect the cable to the connectors on the Ethernet board.

8. Pull in the bottom of the indicator. Re-install the 4 screws (Torque is $1.8N \cdot m \pm 0.5$). Fasten the cap nut of cable gland.

SETUP

Upon installation, the indicator will recognize the Ethernet interface kit and add related items to the menu. Configure the indicator to the desired RS232 and printing parameters (refer to the indicator's instruction manual for assistance).

Ethernet Connection

If DHCP is set to be **ON**, the IP share or router will automatically assign IP Address. If the DHCP is set to be **OFF**, users need to setup the IP Address, Subnet Mask, Gateway, Primary DNS and Secondary DNS. The Interface has a preset IP Address. The Port Number is always **9761** and cannot be changed.

Note: After the setup of the IP Address, Gateway, Preferred DNS and Alternate DNS, please reboot the indicator to enable the setting.

OPERATION

Output

With the interface installed, the indicator will operate in one of the following two ways:

- **Auto Print = OFF, Stable = ON or OFF**
Press **PRINT** will send the display data.
If **Stable** is set to be **ON**, the indicator will wait for a stable reading before sending the stable data.
- **Auto Print = ON, Stable = ON or OFF**
The indicator will automatically send data based on the settings in the menu.
If **Stable** is set to be **ON**, only stable values will be sent.

Data sent from the Interface is in standard ASCII format terminated with a [**CR LF**] (carriage return-line feed). For the output format, please refer to the indicator's instruction manual.

Input

The indicator will respond to various commands sent via the interface adapter. Terminate commands with [**CR LF**].

Commands

For the full list of commands, please refer to the indicator's instruction manual.

CONFIGURATION

The Ethernet option board is connected to the i-DT61XWE indicator through the RS232 interface. The baud rate is 19200.

If the option board is not connected, the submenu will not display. All the setting parameters will be set through the OHAUS ScaleMate software. You can contact an OHAUS authorized dealer for the software.

Ethernet	Reset	no, yes
	IP Address	192.168.1.1-9761
	Alt Print CMD	A~Z, P
	Alt Tare CMD	A~Z, T
	Alt Zero CMD	A~Z, Z
	End	\

For printing template setup, please refer to the following table

Print Setup	Reset	no, yes
	Assignment	Demand , Auto On Stable, Auto On Accept, Interval (seconds), MT-Continuous, OH-Continuous, SICS
	Stable Weight Only [Demand]	Off , On (LFT is forced to be On)
	Mode [Auto On Stable]	Load , Load and Zero
	Time [Interval(seconds)]	1 ~50000
	Checksum [MT-Continuous]	On/ Off
	Select Template	Simple , Custom 1, Custom 2, Custom 3, Custom 4, Custom 5
	End Print3	\

Host Name

The **Host Name** is the name of the Ethernet option board in the network. The default is "OHETHER".

MAC Address

The **Medium Access Control (MAC)** Address of the **Ethernet** option board is not editable.

Port

The Ethernet option board has a TCP server, with the port of 9761 (unmodifiable).

DHCP

If DHCP is set to be **ON**, the IP share or router will automatically assign IP Address. If the DHCP is set to be **OFF**, users need to setup the IP Address, Subnet Mask, Gateway, Primary DNS and Secondary DNS. The Interface has a preset IP Address. The Port Number is always **9761** and cannot be changed.

Note: After the setup of the IP Address, Gateway, Preferred DNS and Alternate DNS, please reboot the indicator to enable the setting.

IP Address

IP is 192.168.1.1 by default.

Subnet Mask

Subnet mask is 255.255.255.0 by default.

Gateway

Gateway is 192.168.1.1 by default.

Primary DNS & Secondary DNS

DNS value is not needed when the system is connected to a LAN, while the DNS value is assigned by the ISP when it is connected to the internet.

Alternative Command

The alternative command of Print, Tare, Zero can be set to 'a' ~ 'z' or 'A' ~ 'Z'.

Note:

1. The submenu will be displayed after the installation of the Ethernet option board.
2. To set the IP address and other parameters, you need to go to the ScaleMate software. To obtain it, contact an authorized OHAUS dealer.

COMPLIANCE

Compliance to the following standards is indicated by the corresponding mark on the product.

Mark	Standard
	<p>This product complies with the applicable harmonized standards of EU Directives 2011/65/EU (RoHS) and 2014/30/EU (EMC). The EU Declaration of Conformity is available online at www.ohaus.com/ce.</p>
	<p>This product complies with the EU Directive 2012/19/EU (WEEE). Please dispose of this product in accordance with local regulations at the collecting point specified for electrical and electronic equipment. For disposal instructions in Europe, refer to www.ohaus.com/weee.</p>
	<p>EN 61326-1</p>

ISED Canada Compliance Statement:

This Class A digital apparatus complies with Canadian ICES-003.

ISO 9001 Registration

The management system governing the production of this product is ISO 9001 certified.

FCC Supplier Declaration of Conformity

Unintentional Radiator per 47CFR Part B
Trade Name: OHAUS CORPORATION
Model or Family identification: Ethernet kit

Issuing Party that Assembled the Product:

Ohaus Instruments (Changzhou) Co., Ltd.
2F, 22 Block, 538 West Hehai Road, Xinbei District, Changzhou
Jiangsu 213022
China
Phone: +86 519 85287270

Responsible Party – U.S. Contact Information:

Ohaus Corporation
7 Campus Drive, Suite 310
Parsippany, NJ 07054
United States
Phone: +1 973 377 9000
Web: www.ohaus.com

FCC Compliance Statement:

Note: This equipment has been tested and found to comply with the limits for a Class A digital device, pursuant to part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference when the equipment is operated in a commercial environment. This equipment generates, uses, and can radiate radio frequency energy and, if not installed and used in accordance with the instruction manual, may cause harmful interference to radio communications. Operation of this equipment in a residential area is likely to cause harmful interference in which case the user will be required to correct the interference at his own expense.

Changes or modifications not expressly approved by the party responsible for compliance could void the user's authority to operate the equipment.

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P/N 30572926 A © 2020 Ohaus Corporation, all rights reserved / todos los derechos reservados/ tous droits réservés/ Alle Rechte vorbehalten / Tutti i diritti riservati / 版权所有
OHAUS Corporation