Navigator™
Ethernet Interface
INSTRUCTION MANUAL

Interfase Ethernet
Manual de instrucciones

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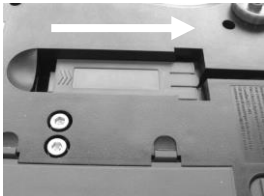
INTRODUCTION



This Ethernet interface kit is for use with the Ohaus Navigator Series Scales only. This accessory has been tested and complies with the approvals listed in the Navigator Instruction Manual.

INTERFACE INSTALLATION

1. Disconnect the existing Power Adapter, it will be replaced.
2. Remove the blank interface cover by sliding the cover back.

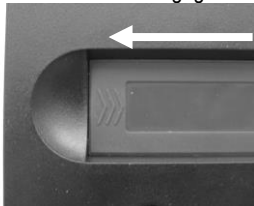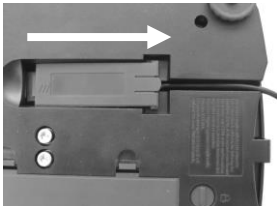3. Install the Interface Module. Slide the cover forward to engage the connections.

4. Connect an Ethernet Cable to the RJ45 connector.

SETUP

Upon installation, the scale will recognize the Ethernet Interface Kit and add the following items to the menu structure.

CAL	SETUP	UNITS	PRINT	Ethrnt	END
			Stable	On-Off	
			On, Off	On, Off	
			A.Print	IP Adr	
			On.stbl, On.Acc*,	End	
			Cont, 5sec, 15sec,		
			30sec, 60sec, off		
			End		

*Note: Print and Ethernet menu selections may vary between the different balance series.

Setup Definitions

PRINT / Stable - On	Transmits stable data only
PRINT / A.Print – On.Stbl	Automatically outputs data upon stability
PRINT / A.Print – On.Acc	Automatically outputs data in Accept Range
PRINT / A.Print – Cont	Repeatedly sends data as fast as possible
PRINT / A.Print – (xx) SEC	Sends data every (xx) seconds
PRINT / A.Print – OFF	Sends data when PRINT is pressed or P command
E.t.h.r.n.t / E.t.h.r.n.t - OFF	Turns Interface off, saving battery power
E.t.h.r.n.t / Ip Adr	Displays the scales IP Address

Ethernet Connection

The Interface has a preset IP Address which can be displayed by entering **Ip Adr** in the menu. Use the “No” and “Back” keys to show the entire number on the display. The Port Number is **9761**.

OPERATIONOutput

With the interface installed the balance will operate in one of the following three ways:

- Ethernet = on, Auto Print = off, Stable = on or off
Pressing PRINT will send the display data. If Stable is set to On, the balance will wait for a stable reading before sending data.
- Ethernet = on, Auto Print = on, Stable = on or off
The balance will automatically send data based on the settings in the Menu. If Stable is set to On, only stable will be sent.
- Ethernet = off
During battery operation, turning the Ethernet Interface off saves significant power.

Data sent is in standard ASCII format terminated with a [CRLF] (carriage return-line feed).

[weight]	10 characters	(right justified)
[space]	1 character	
[unit]	5 characters max	(left justified)
[space]	1 character	
[stability indicator]	1 character	“?” when unstable, blank when stable
[CR]	1 character	
[LF]	1 character	

Output examples: (Note: * and _ indicate spaces, CR/LF not shown)

****192.21_g	- Manual, Continuous or On-Stable printing
*****0.01_g_?	unstable reading
*****0.01_g__00:00:00	- Interval Printing (example:15sec interval)
****176.30_g_?_00:00:15	unstable reading
****192.08_g__00:00:30	
****192.21_g	
****207.80_g_TOTAL	- Accumulate Mode (Manual printing only)

Input

The scale will respond to various commands sent via the interface. The commands must be terminated with a [CR] or [CRLF].

OP	Turns Auto-Print off
SLP	Auto-Print on, prints on stability (non-zero value only)
SLZP	Auto-Print on, prints on stability (zero and non-zero values)
CP	Continuous Auto-Print
(x)P	Auto-Print on 1 to 3600 second intervals (x = 1 to 3600)
PM	Sends current mode
M	Advance to next enabled mode
PU	Sends current unit
U	Advance to next enabled unit
T	Tares balance, same as pressing TARE
Z	Zeros balance, same as pressing ZERO
PV	Sends software version
P	Same as pressing PRINT
IP	Immediately sends stable or unstable value
SP	Sends stable data only

Auto-Print Operation

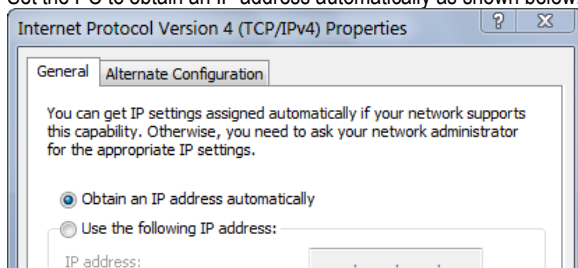
Once Auto-Print is activated in the menu, the balance will send data as required. To temporarily stop Auto-Printing, press the PRINT key. If there is data in the print buffer the printer will finish printing this data. A second press will resume Auto-Printing.

(Note: In some modes, stopping or resuming Auto-Print must be done when the balance is not reading "0".)

ETHERNET CONFIGURATION

To set the Ethernet parameters per web.

1. Make sure the Ethernet Interface Kit is installed properly and network cable is connected to a computer (PC).
2. Set the PC to obtain an IP address automatically as shown below.

- Open the IE Browser and enter the default IP address of the Ethernet Interface Kit

"169.254.1.1" () , the browser will show below page:

- Click the "Network Configuration" on the left side, and input the user & password. Then click "OK" to proceed.
- Input the Ethernet parameters in the following page, click "Save Config" to save the changes.

MAC Address:	00:E0:7C:00:00:00
Host Name:	OHAUETH
<input checked="" type="checkbox"/> Enable DHCP	
IP Address:	169.254.1.1
Gateway:	169.254.1.1
Subnet Mask:	255.255.0.0
Primary DNS:	169.254.1.1
Secondary DNS:	0.0.0.0
<input type="button" value="Save Config"/>	

Note:

- To use DHCP, enable the DHCP: **Enable DHCP**
- To use static IP address, disable the DHCP: **Enable DHCP** and then set the Ethernet configuration parameters.

ACCESSORIES

For a complete listing of Ohaus printers and other accessories, contact Ohaus Corporation or visit www.Ohaus.com

COMPLIANCE

This accessory has been tested and complies with the approvals listed in the Navigator Instruction Manual.

Disposal

In conformance with the European Directive 2012/19/EU on Waste Electrical and Electronic Equipment (WEEE) this device may not be disposed of in domestic waste. This also applies to countries outside the EU, per their specific requirements.

Please dispose of this product in accordance with local regulations at the collecting point specified for electrical and electronic equipment. If you have any questions, please contact the responsible authority or the distributor from which you purchased this device.

Should this device be passed on to other parties (for private or professional use), the content of this regulation must also be related.

For disposal instructions in Europe, refer to www.ohaus.com/weee.

Thank you for your contribution to environmental protection.

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P/N 83032209 B © 2018 Ohaus Corporation, all rights reserved / todos los derechos reservados / tous droits réservés / Alle Rechte vorbehalten / tutti i diritti riservati / все права защищены.